

# MAX300-CAT Startup Guide

824285 Revision 1  
February 2020

## I. Connecting the Hardware

1. Connect the serial cable provided to the MAX300-CAT Mass Spectrometer Electronics



2. Connect other end of the serial cable to the USB to serial converter provided and connect the USB to the computer




3. If a stream selector option was purchased, connect the gray serial cable from the MAX300 CAT stream selector to the serial port on the computer



4. Power ON the Pumping station




5. Press the START button (F1) on the pump.

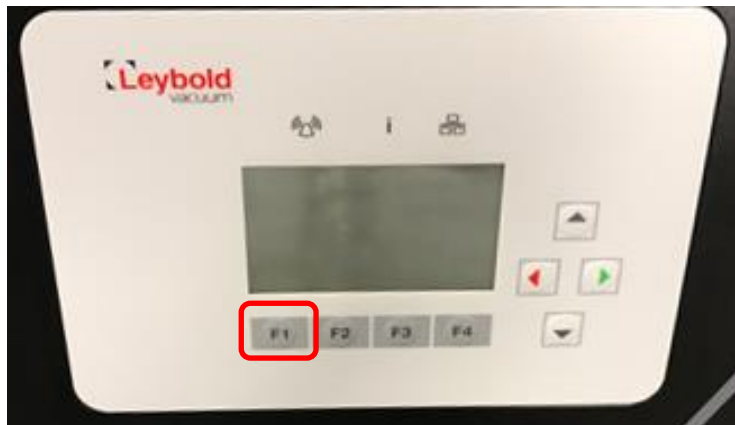


## II. Install the USB to RS-232 Serial Converter Driver.

For Windows 7: (See next section for Windows10 instructions)


1. Power on the Computer
2. Right-click "My Computer" and select "Manage"
3. Click "Device Manager"

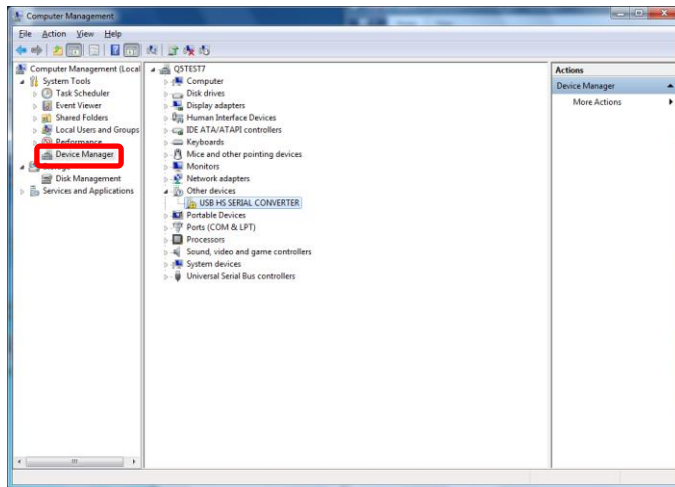


4. Right-click "USB HS SERIAL CONVERTER" and select "Update Driver Software"
5. Choose "Browse my computer for driver software"

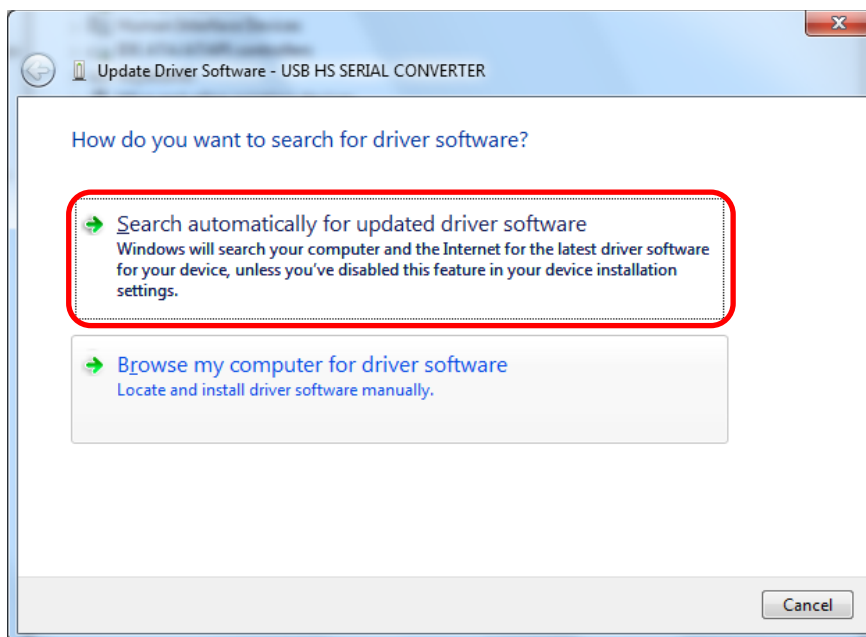


6. Click the "Browse" button

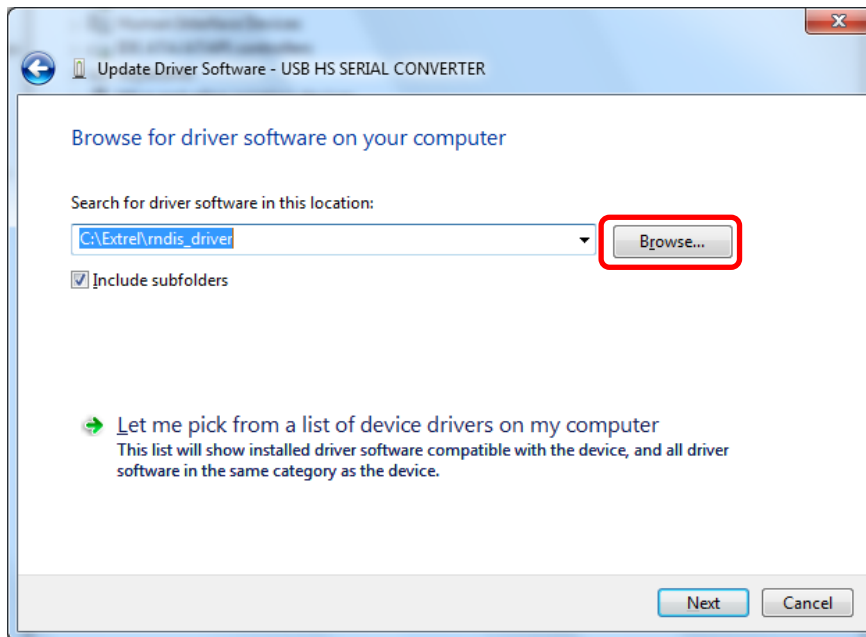


7. Choose the folder "i386" inside the "USB to serial Converter Driver" folder on the USB disk provided

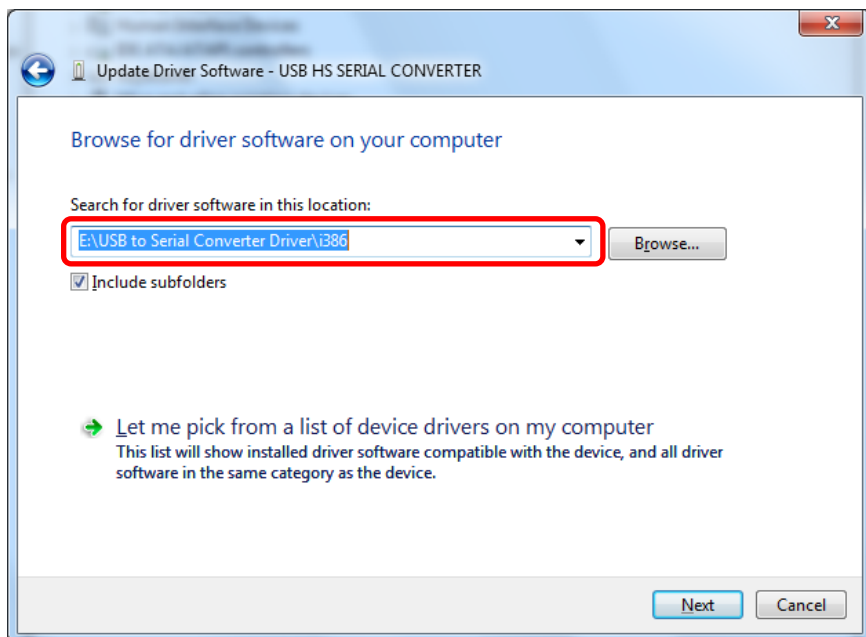


8. Check the box labeled "Always trust software from..." and click the "Install" button

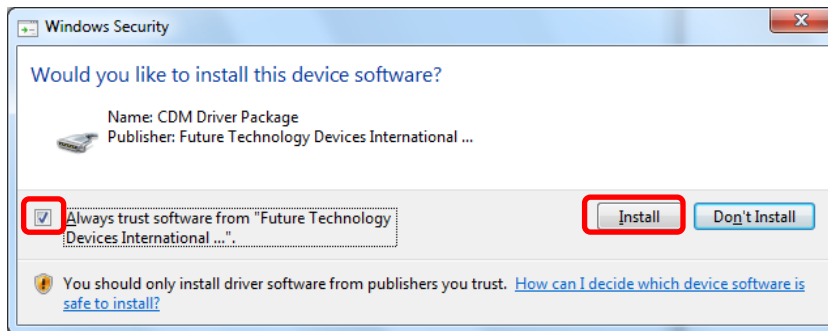


9. Allow the software to complete the installation

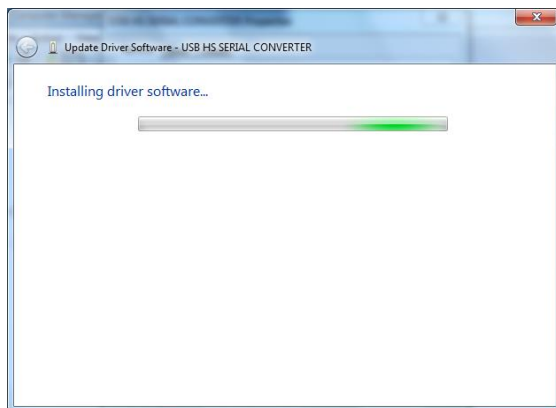

10. Click the "Close" button

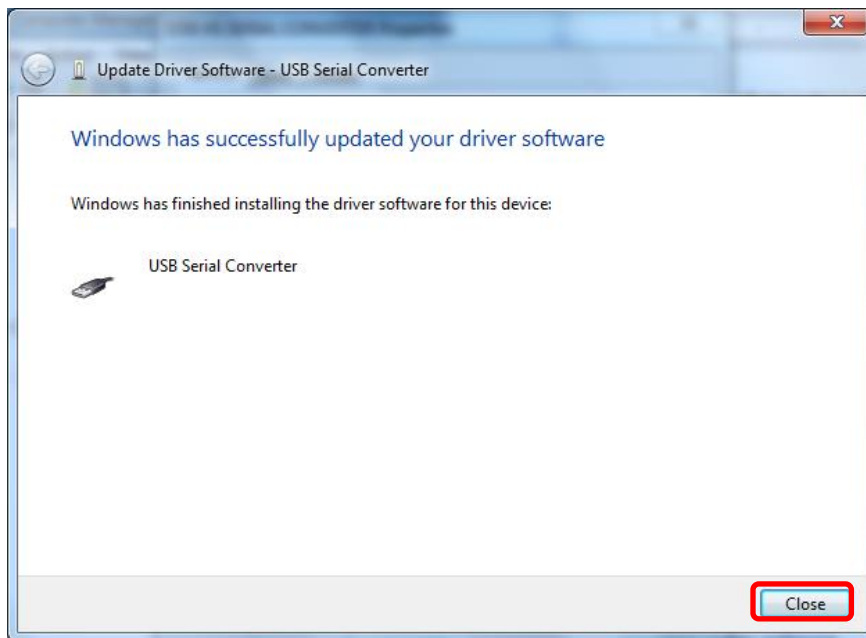


**\*\*NOTE** which COM port is labeled "USB Serial Port"

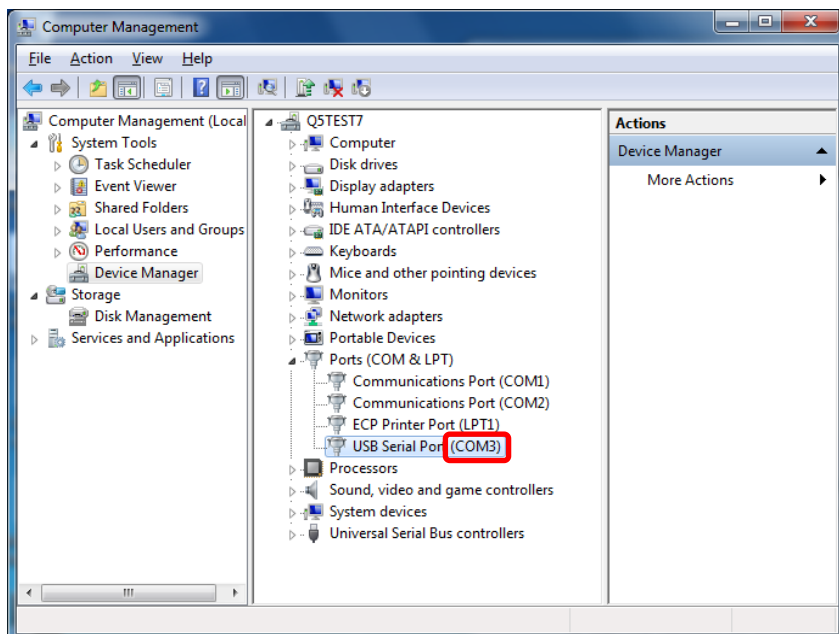


## USB to Serial Converter driver installation for Windows10

**Note: The Computer should have a valid Internet connection to complete this step.**



1. Power on the Computer
2. Right-click "This PC" and select "Manage"
3. Click "Device Manager"

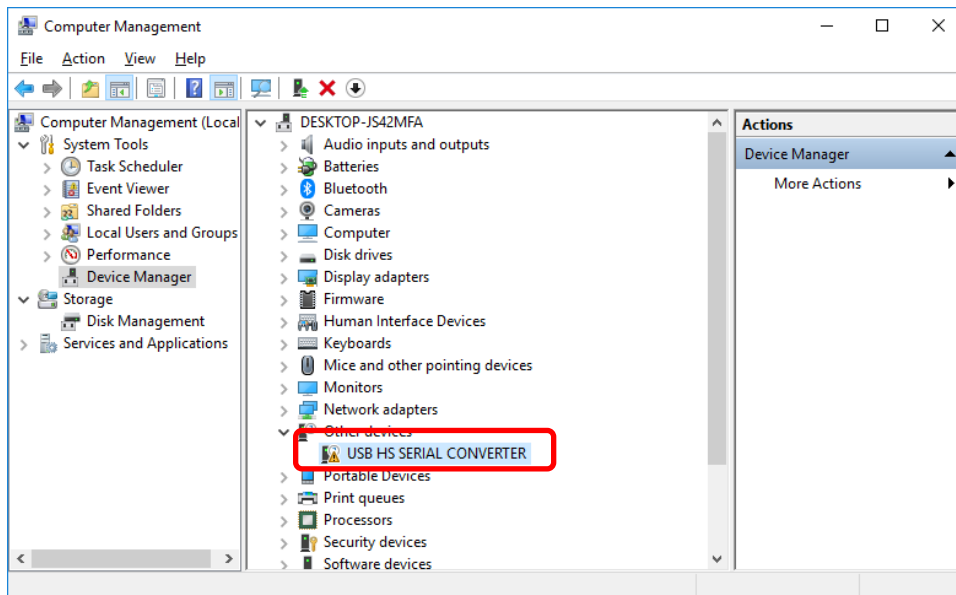


4. Right on "USB HS SERIAL CONVERTER" and select "Update Driver"

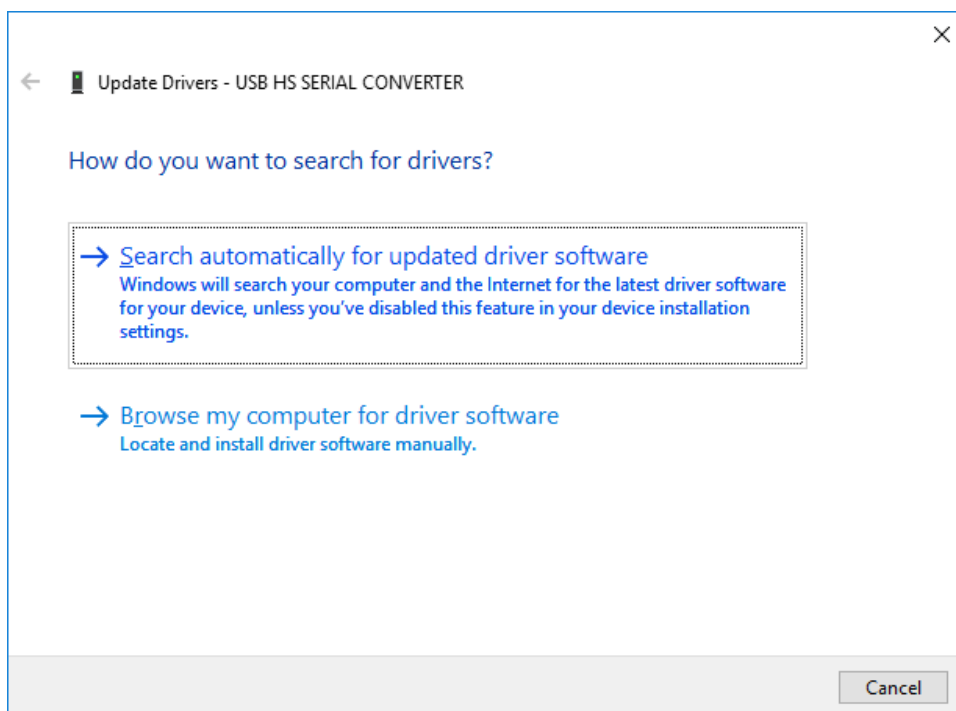



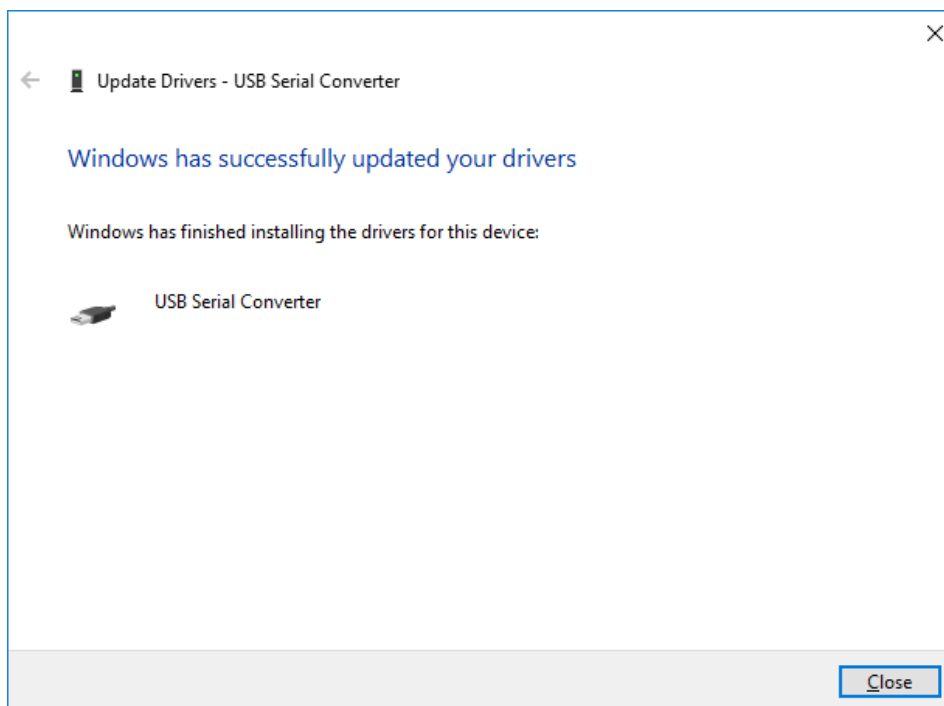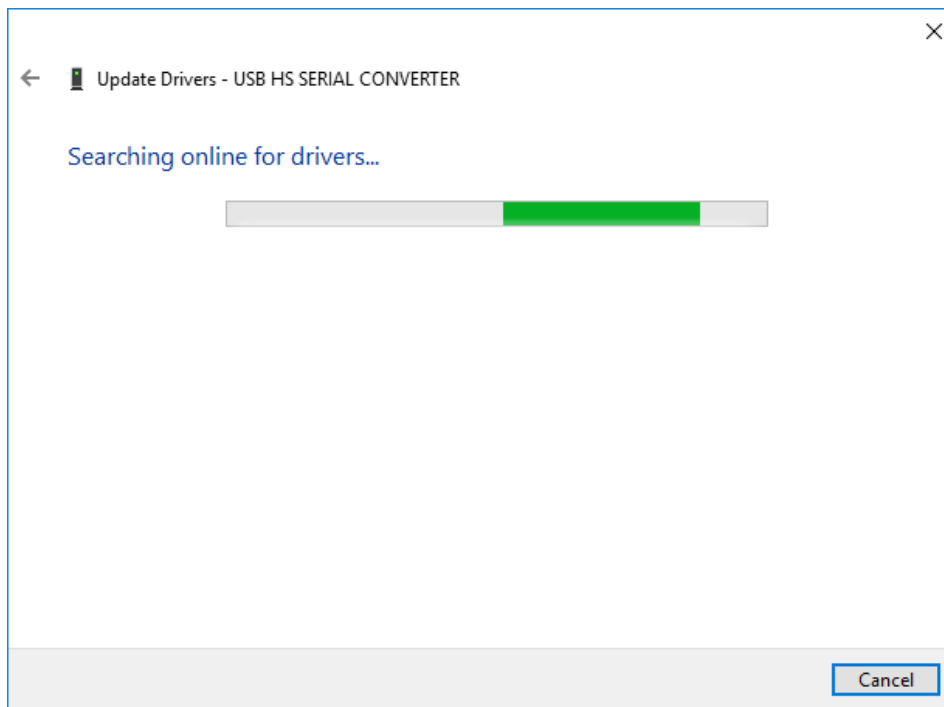


**\*\*Note down which COM port is labeled "USB Serial Port". This port name must be specified in the instrument configuration file as described in a later step.**

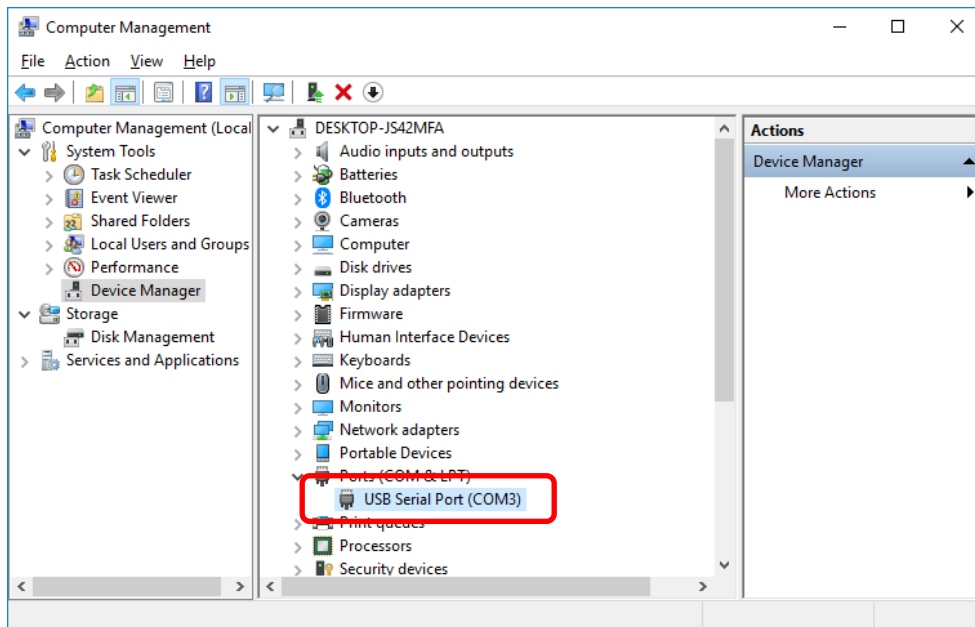

### III. PC Setup

Make sure the Power options are set in the Windows so that the PC would never goes to sleep mode. If the PC goes to sleep while collecting scan data, system will be stalled and collection stops.

You can access the Power options window from Control Panel – System and Security – Power Options. Set the options to “Never” sleep as shown below:

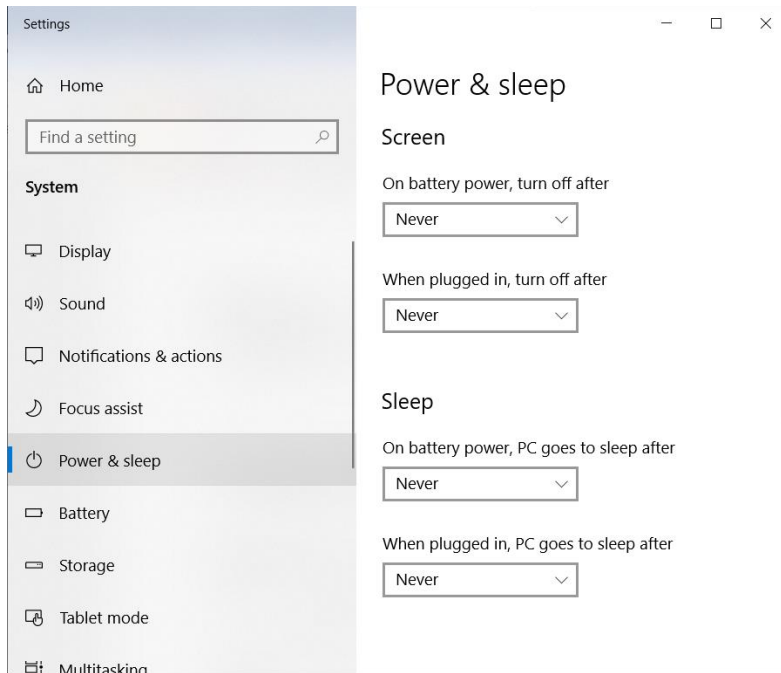


### IV. Software Installation and Setup

#### Install the Questor5 Software

8. Locate the “pre\_install.bat” file in the “Questor5” folder on the USB disk provided

Right-click & select "Run as Administrator" option


9. When the installer finishes, press any key to exit.

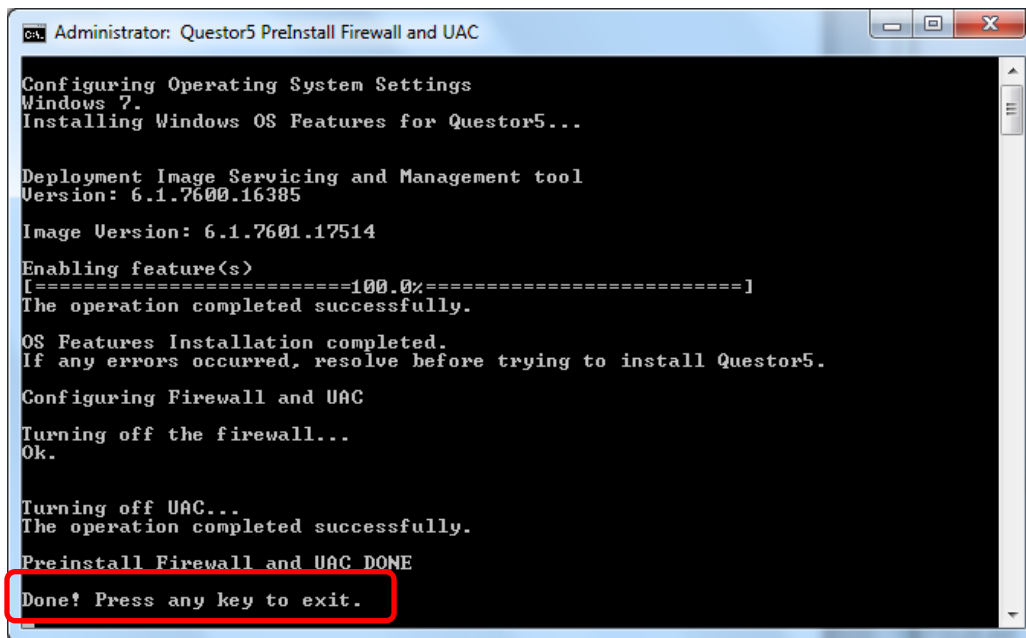


10. Locate the “setup.exe” file in the “Questor5 folder on the USB disk provided

Right-click & select "Run as Administrator" option


11. When prompted, select the “Install” button

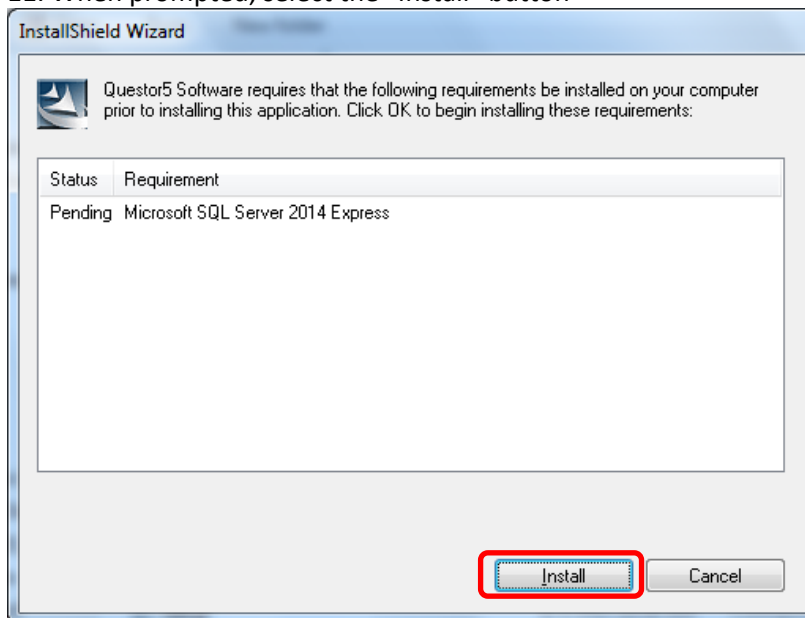


12. Click “Next”

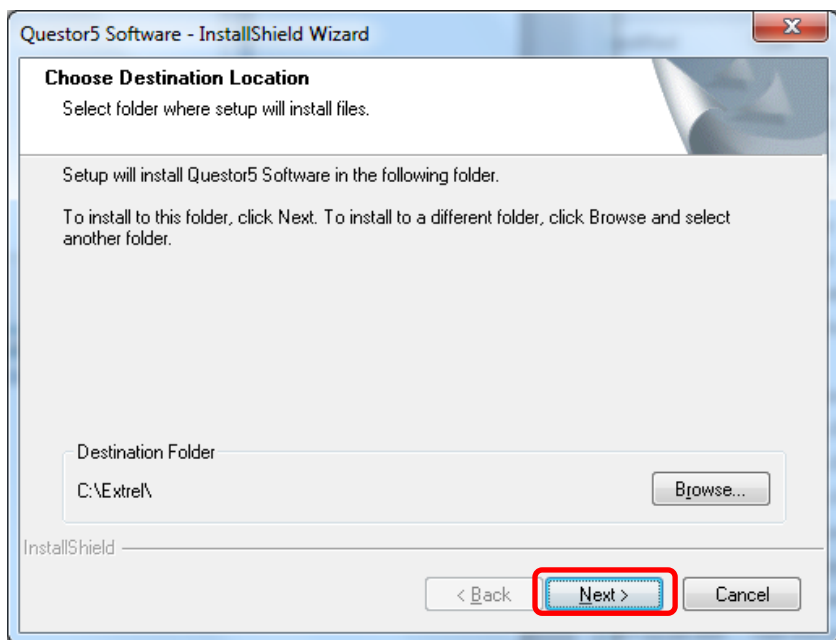


13. Choose “MAX300 CAT / RGA”

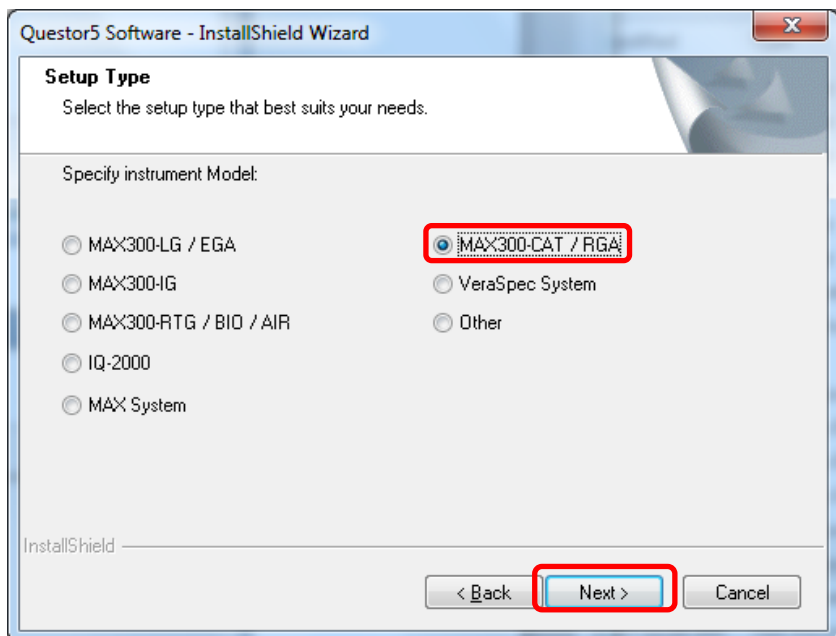


14. Select the sample selector option purchased with the system (if no sample selector option was purchased, choose “No Sample Selector installed”)

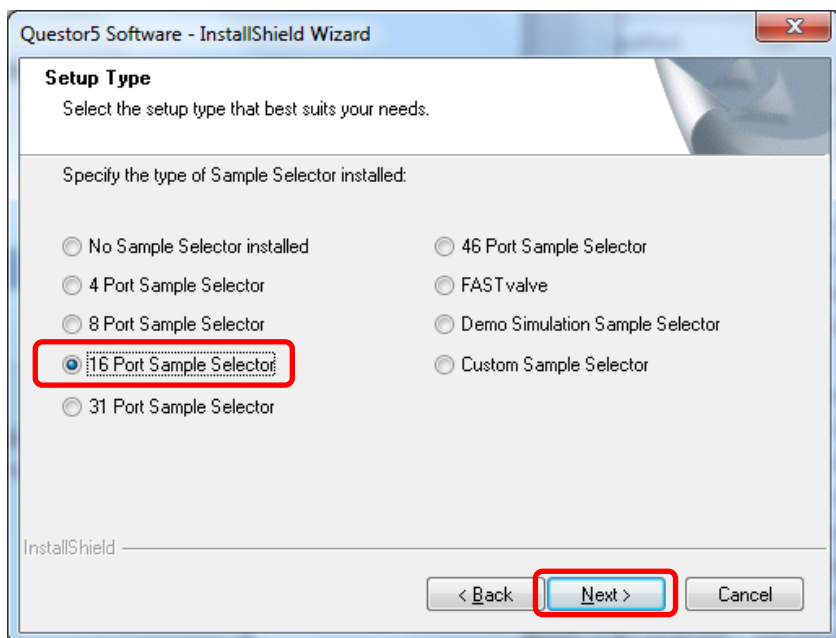


15. Click "Next"

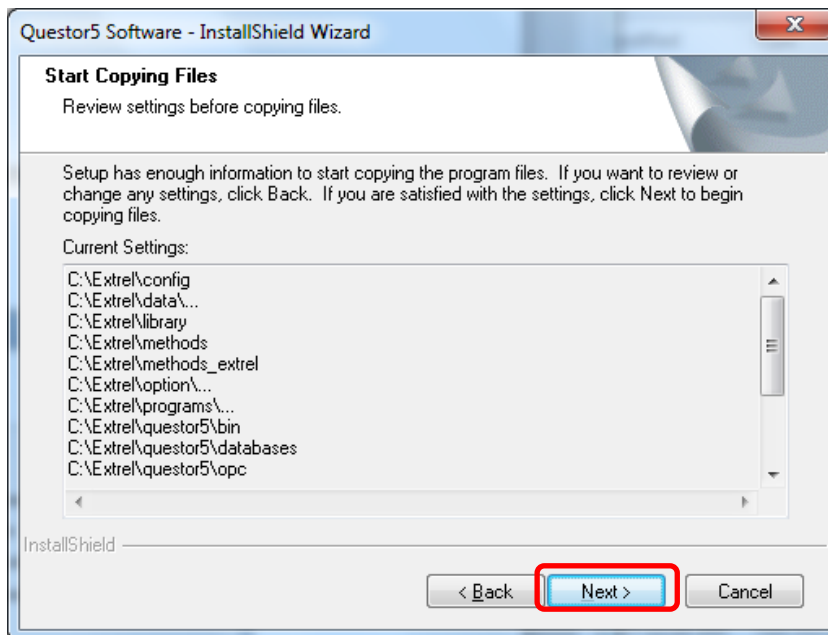

16. Wait for the software to complete its installation

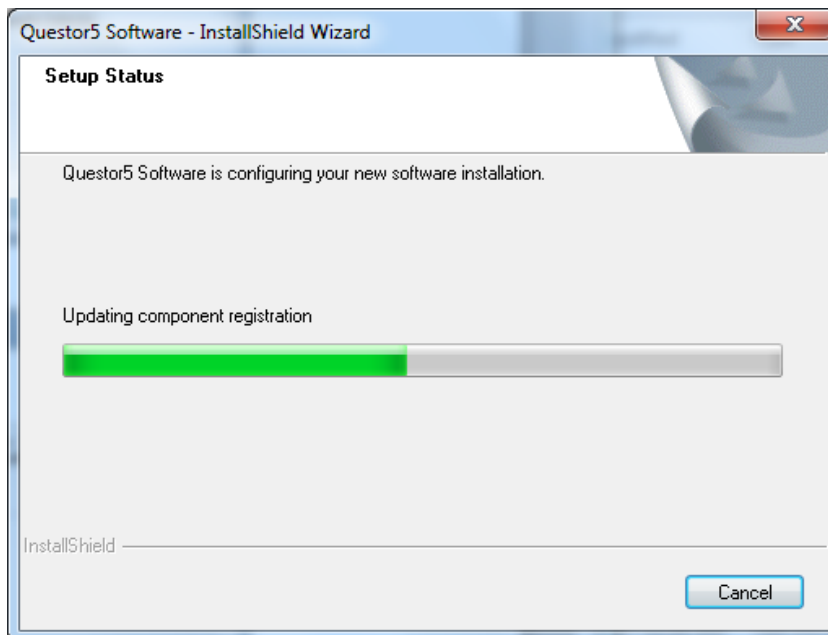


17. Click "Finish" to complete the installation and restart the computer

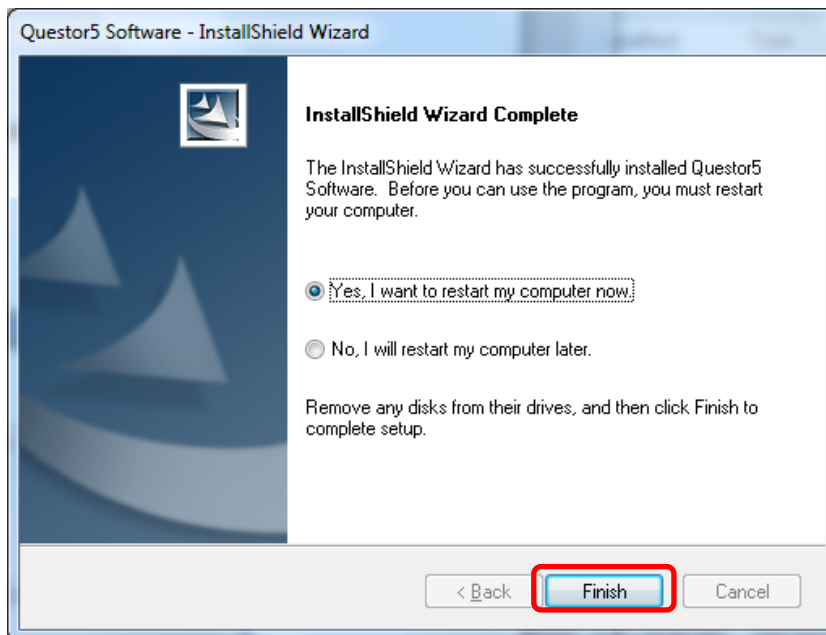

18. Locate the "post\_install.bat" file in the "Questor5" folder on the USB disk provided

Right-click & select "Run as Administrator" option

19. When the installer finishes, press any key to exit

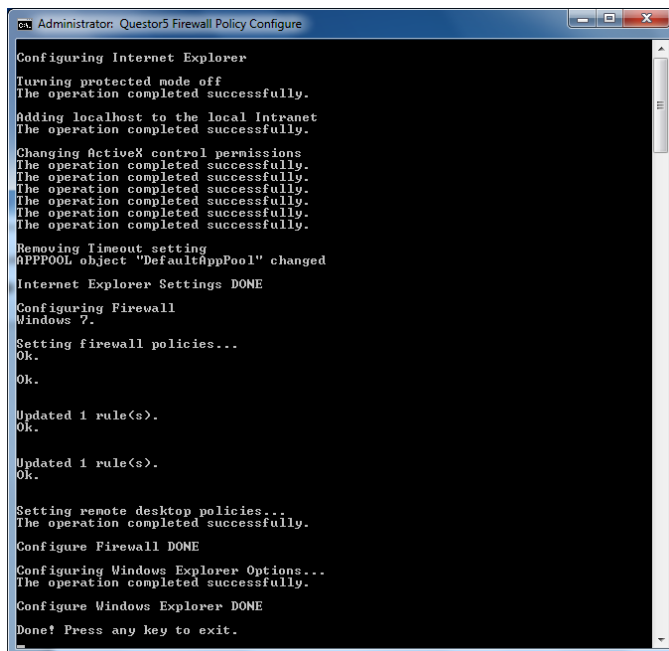


20. Reboot the computer



## C. Software Setup

1. Copy the Configuration File for the instrument "sn####\_factory\_cal.cfg" from the "Customer Files" folder on the USB disk to the C:\Extrel\config folder.
2. Open the above file in notepad and modify the "Port" attribute to the same port name noted during the USB serial converter driver installation.
3. Change the path in the "Filename" attribute to "C:\Extrel\Config\" as shown below:

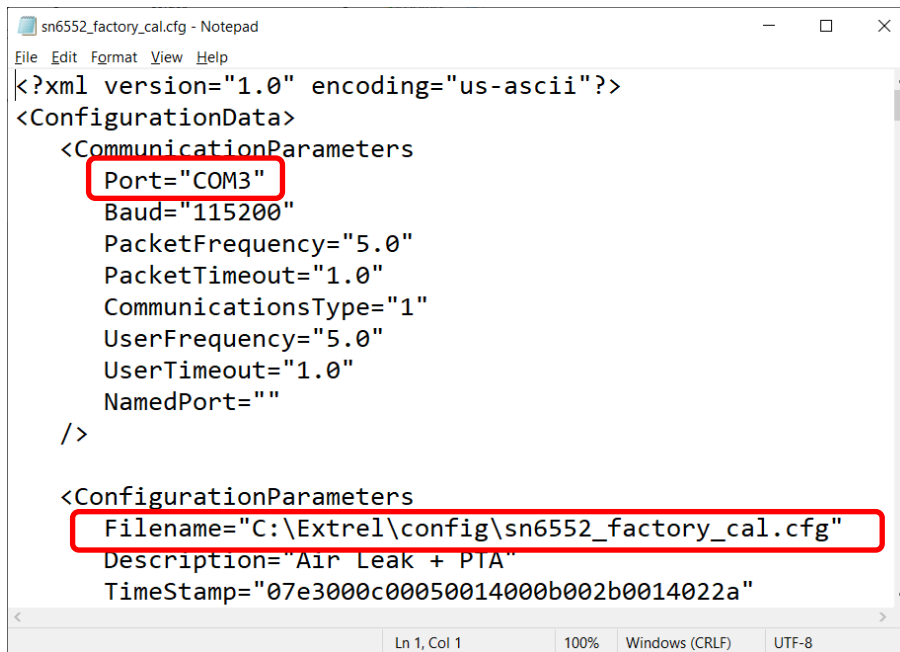

```
sn6552_factory_cal.cfg - Notepad
File Edit Format View Help
<?xml version="1.0" encoding="us-ascii"?>
<ConfigurationData>
  <CommunicationParameters
    Port="COM3"
    Baud="115200"
    PacketFrequency="5.0"
    PacketTimeout="1.0"
    CommunicationsType="1"
    UserFrequency="5.0"
    UserTimeout="1.0"
    NamedPort=""
  />
  <ConfigurationParameters
    Filename="C:\Extrel\config\sn6552_factory_cal.cfg"
    Description="Air Leak + PIA"
    TimeStamp="07e3000c00050014000b002b0014022a"
  />
/>
Ln 1, Col 1    100%    Windows (CRLF)    UTF-8
```

4. In the same file, make sure the "ScanSpeed" is set to 144, "SamplesPerAMU" to 20 and "FilamentForceOffStartup" to 1 as highlighted below:

```
sn6552_factory_cal.cfg - Notepad
File Edit Format View Help

<ScanParameters
  ScanSpeed="144"
  LowMass="1"
  HighMass="310"
  SamplesPerAMU="20"
/>

<OperatingParameters
  Mode="Mass sweep"
  Focus1="-22"
  Focus2="-68"
  ElectronEnergy="70"
  FilamentEmission="2.0"
  AutoZero="Off"
  ScanMode="Sweep"
  Filament="1"
  PressureUnits="Torr"
  EnableElectronMultiplier="0"
  MultiplierVoltage="1020"
  FilamentForceOffStartup="1"
/>

Ln 15, Col 21 100% Windows (CRLF) UTF-8
```


5. Keep a note on the highlighted values, this may need to apply while tuning the instrument.

```
sn6552_factory_cal.cfg - Notepad
File Edit Format View Help

<CalibrationParameters
  LowCalMass="1"
  LowCalResolution="610"
  LowCalPosition="0.42"
  LowCalIonEnergy="5.0"
  HighCalMass="300"
  HighCalResolution="2345"
  HighCalPosition="0.50"
  HighCalIonEnergy="5.0"
  TotalAmpOffset="2000"
  PartialAmpOffset="2200"
  TotalIntegratingCap="10.00"
  PartialIntegratingCap="3.00"
  RFSettleTime="50"
  SWSettleTime="10"
  Pirani1ATM="1.86300"
  PiraniZero="0.28700"
  PiraniAutoRecalibrate="0"
  PartialSensitivity="6.00e-004"
  TotalSensitivity="1.00e+001"
  debug="0"
/>

Ln 1, Col 1 100% Windows (CRLF) UTF-8
```

6. Open the file ExtrelRga3.ini from “C:\Extrel\Programs” folder in notepad and modify the attribute “rga\_cfg” to point to the cfg file referenced in step 1.

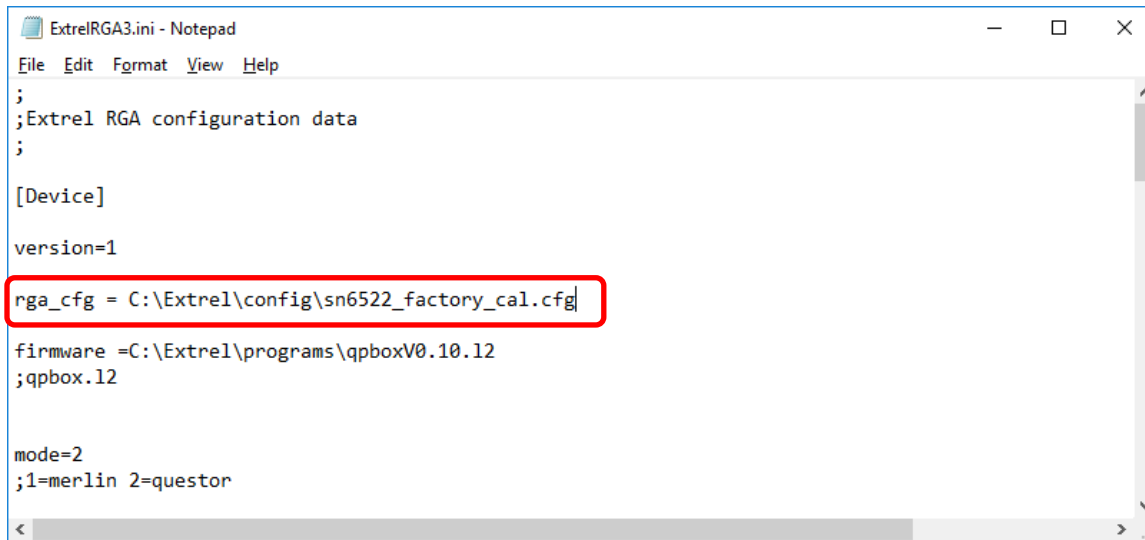A screenshot of a Notepad window titled "ExtrelRGA3.ini - Notepad". The window shows the following text: 


```
;
;Extrel RGA configuration data
;

[Device]
version=1
rga_cfg = C:\Extrel\config\sn6522_factory_cal.cfg
firmware =C:\Extrel\programs\qpboxV0.10.12
;qpbox.12

mode=2
;1=merlin 2=questor
```

 The line "rga\_cfg = C:\Extrel\config\sn6522\_factory\_cal.cfg" is highlighted with a red rectangular box.

7. In the same file, change the “rga\_last\_mass” to 100 or 300 based on the mass range of your system as shown below:

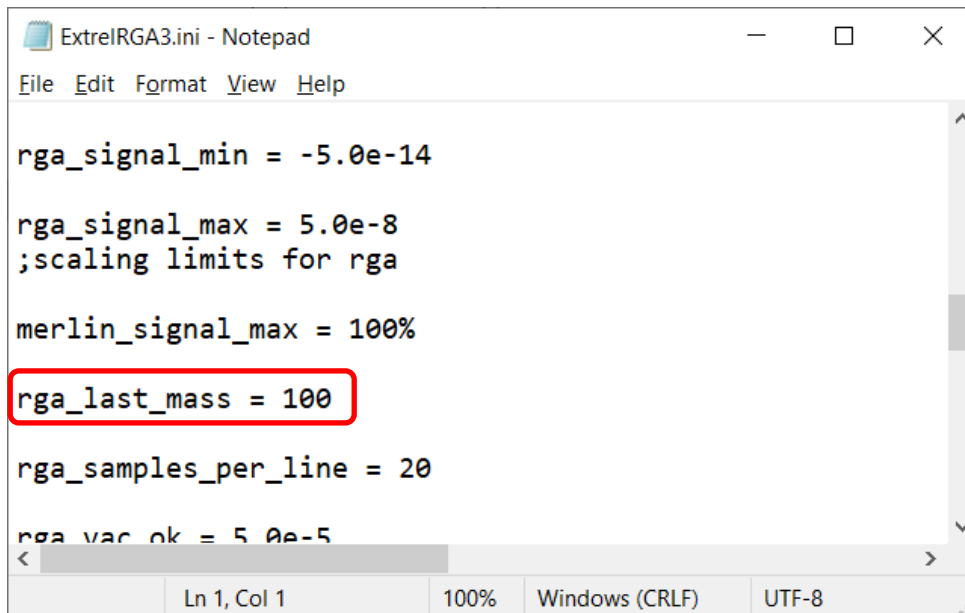A screenshot of a Notepad window titled "ExtrelRGA3.ini - Notepad". The window shows the following text: 


```
rga_signal_min = -5.0e-14
rga_signal_max = 5.0e-8
;scaling limits for rga
merlin_signal_max = 100%
rga_last_mass = 100
rga_samples_per_line = 20
rga_vac_ok = 5.0e-5
```

 The line "rga\_last\_mass = 100" is highlighted with a red rectangular box. The status bar at the bottom of the window displays "Ln 1, Col 1", "100%", "Windows (CRLF)", and "UTF-8".


8. Select the Questor5 Software from the Desktop.




9. Install ActiveX controls by clicking the "Yes" button

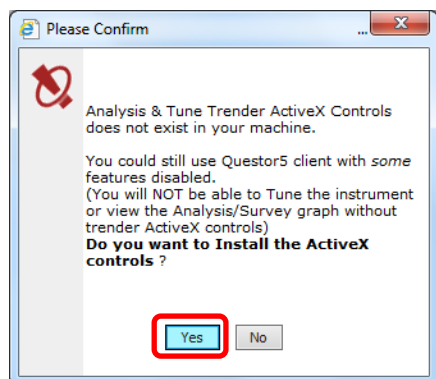


10. Choose "Click here" to install ActiveX controls

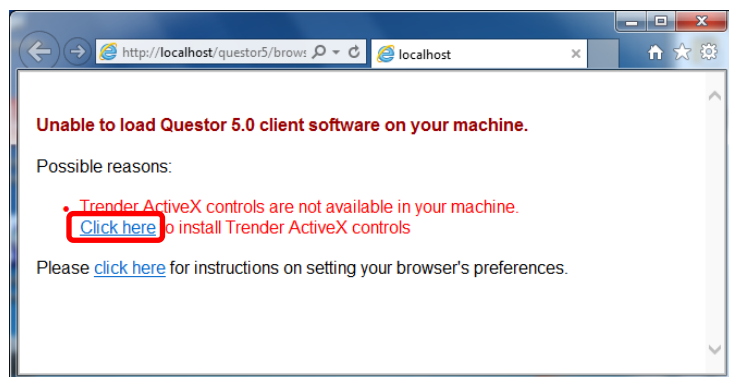


11. Click "Boot Server"

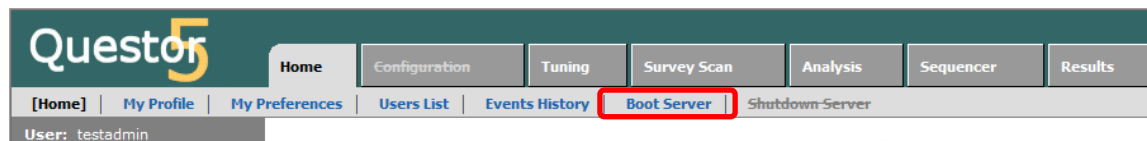


12. Ensure the instrument is online by clicking "Instrument Status"

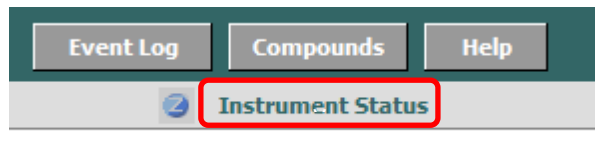


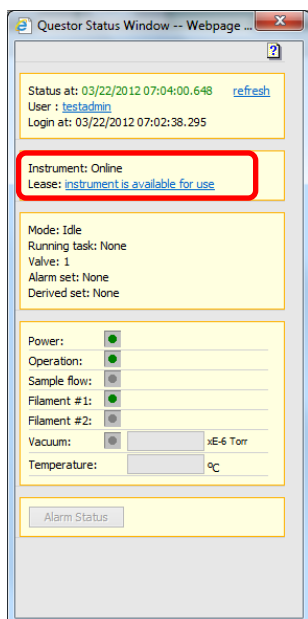


13. Select the "Tuning" tab, and click "Create New"

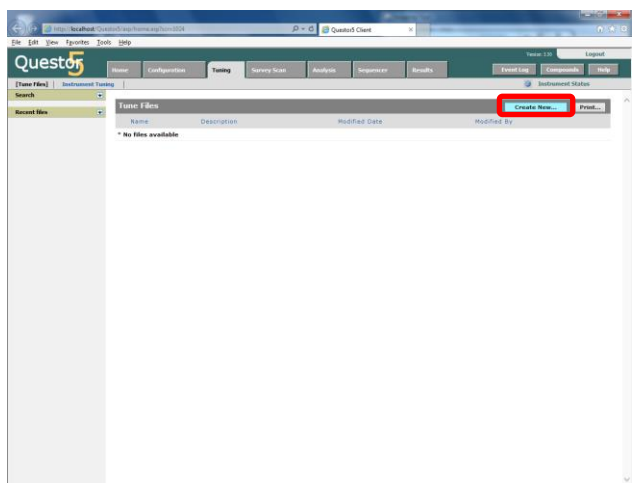


14. Name the Tune File and click "OK"

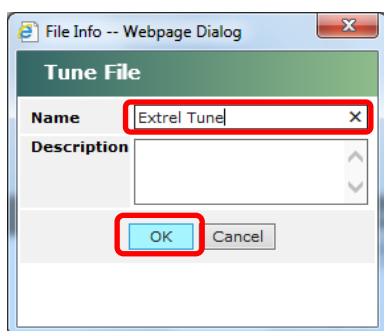

15. Select a mass to start tuning and click "OK"

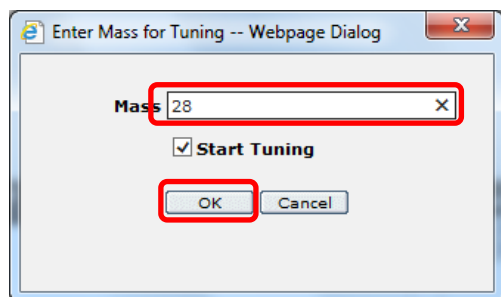


16. Change the tuning start and end masses from 27.5 and 28.5 to 1 and 50, respectively and click "Apply" button. If this is the first time that the instrument has been turned on allow it to run in this mode for at least **3 hours** before attempting to fine-tune the peaks.

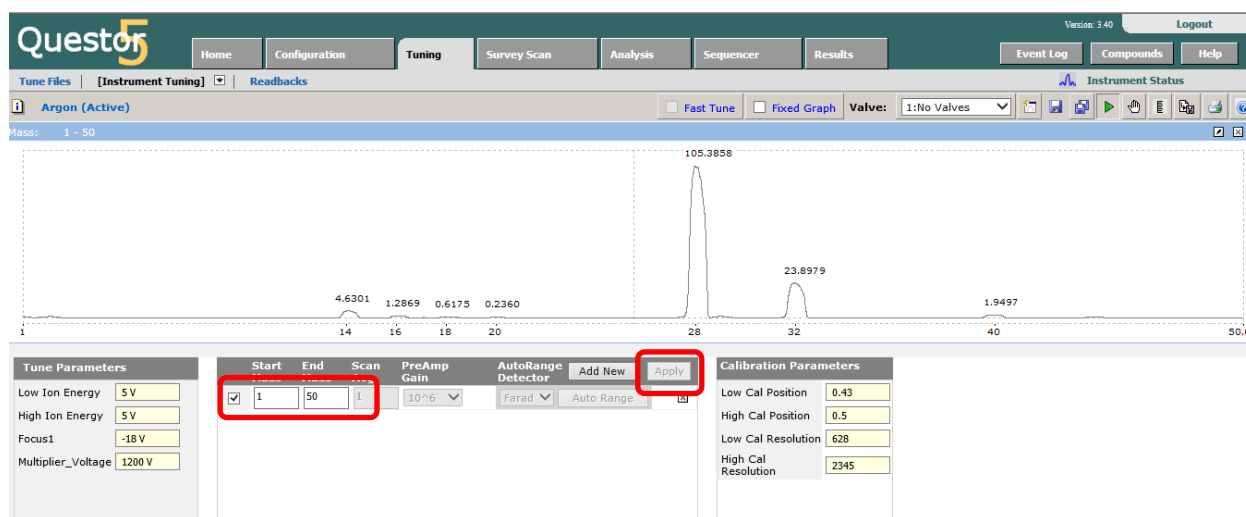


17. In order to accurately tune the peak position and width, create multiple mass ranges 2 amu wide by clicking the "Add New" button. For example, if the instrument is running in air these mass ranges can be useful: 13 – 15 amu, 27 – 29 amu, 31 – 33 amu and 39 – 41 amu. Display these mass ranges and the corresponding peaks side-by-side by clicking the check box to the left of the mass range as shown in the Figure below.

First the resolution must be adjusted so that the peaks are approximately 1 amu wide at their base. To check the width place the mouse to the left and right of each peak and the mass position will be shown in the upper right corner of the corresponding window. The peak width can be adjusted by changing the "Low Cal Resolution" and "High Cal Resolution" parameters which, as the names say, affect more the low and high mass peaks, respectively.


After the resolution has been tuned the peak positions can be adjusted by changing the "Low Cal Position" and "High Cal Position" parameters.

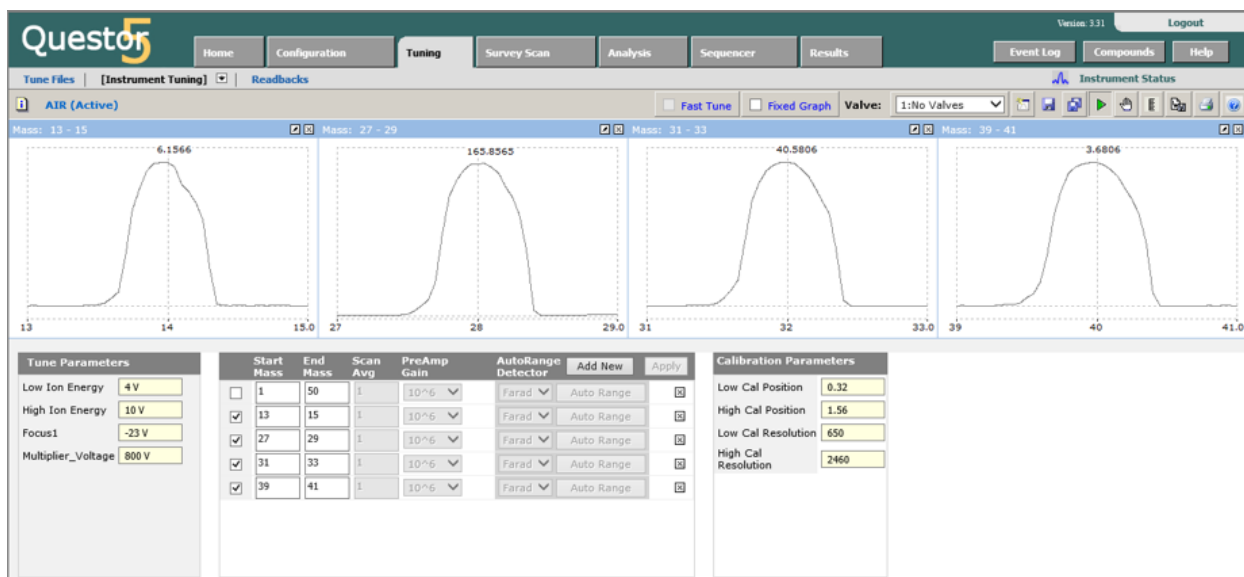

18. Create an analysis file by selecting the “Analysis” tab and clicking “Create New”

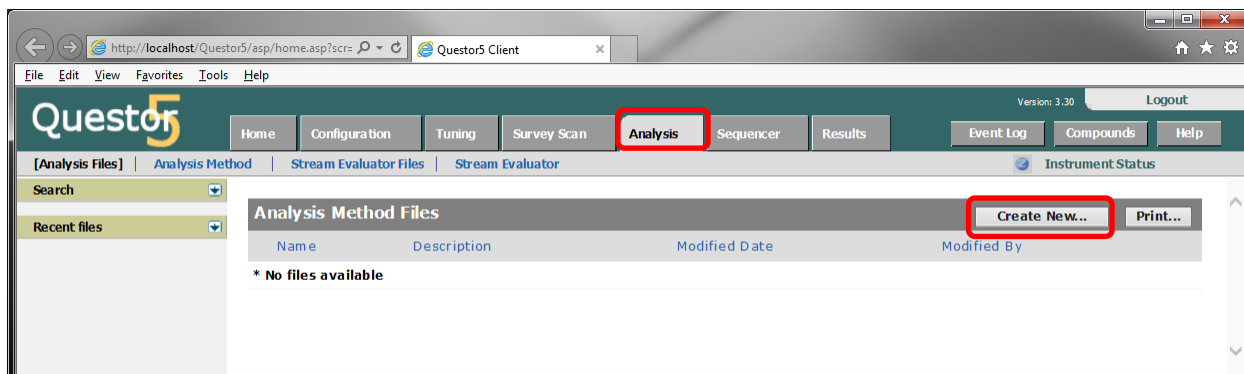


19. Name the Analysis File and click “OK”

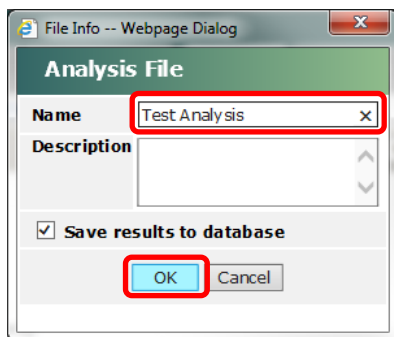


20. Select each of the components in the analysis sample or experiment, one at a time, from the Chemicals List and click the “Select” button. A message will appear near the top of the window saying that the component has been added to the fragment matrix. Once all components have been selected,

click the “Close” button.

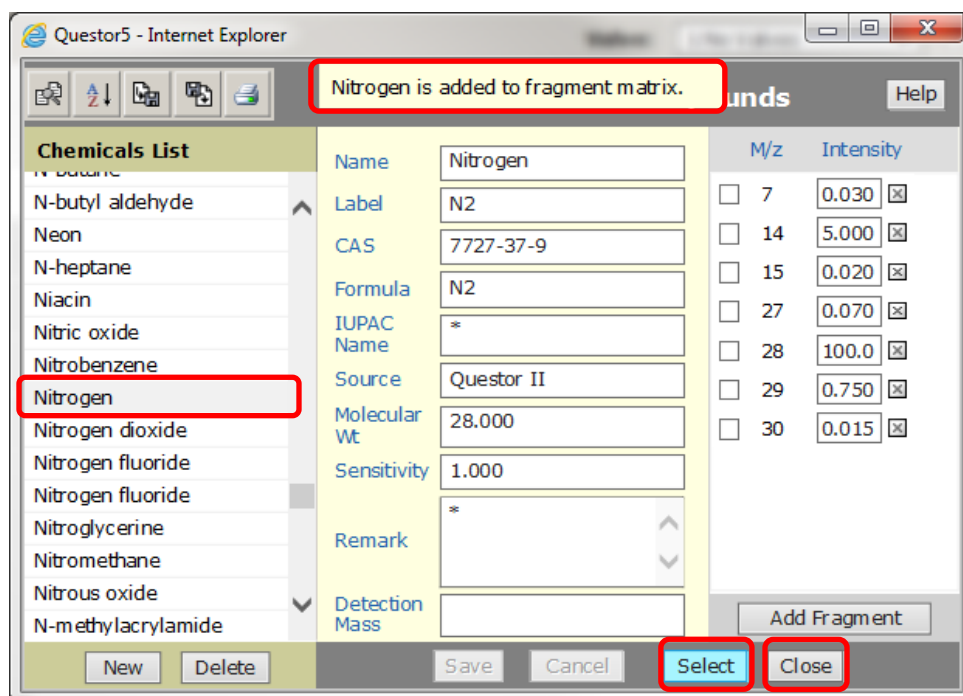

21. Click the “Save” button to save the analysis

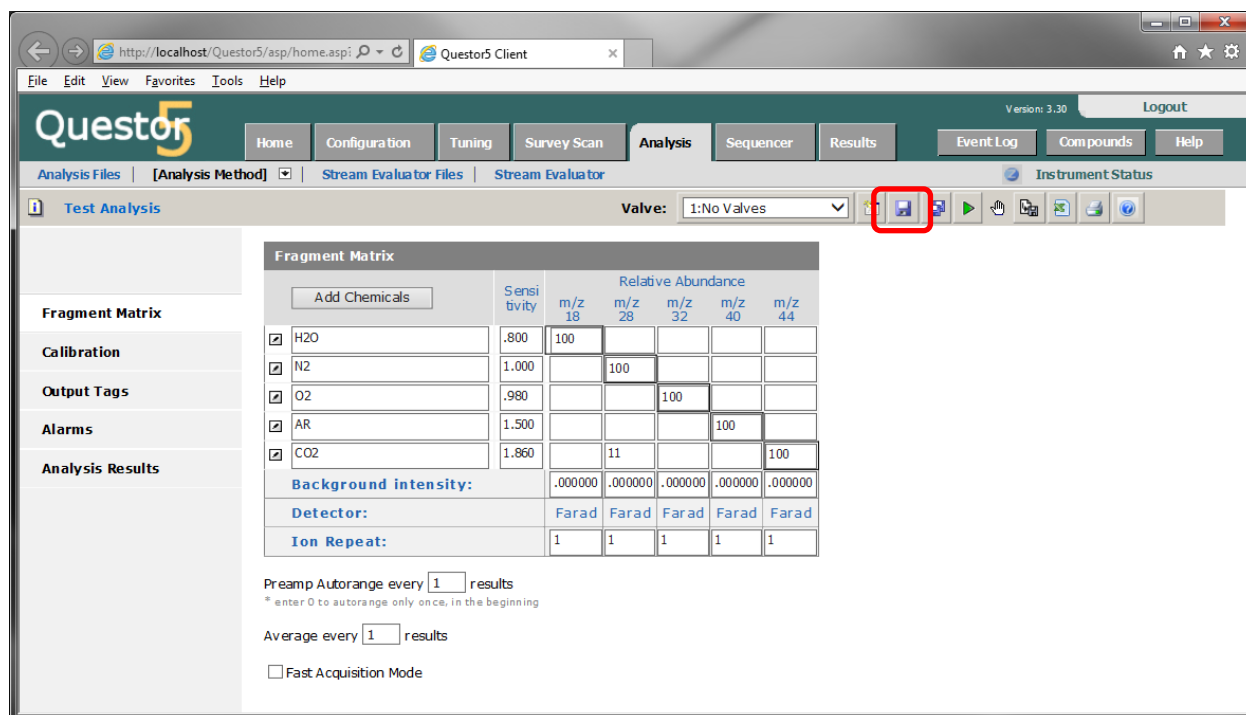

22. If applicable, choose the appropriate valve position from the “Valve” dropdown menu and click the “Start Analysis” button



Questor5 Client

Version: 3.30 Logout

Home Configuration Tuning Survey Scan **Analysis** Sequencer Results Event Log Compounds Help

Analysis Files [Analysis Method] Stream Evaluator Files Stream Evaluator Instrument Status

Test Analysis Valve: 1:No Valves

**Fragment Matrix**

Add Chemicals

	Sensitivity	Relative Abundance				
		m/z 18	m/z 28	m/z 32	m/z 40	m/z 44
<input checked="" type="checkbox"/> H2O	.800	100				
<input checked="" type="checkbox"/> N2	1.000		100			
<input checked="" type="checkbox"/> O2	.980			100		
<input checked="" type="checkbox"/> AR	1.500				100	
<input checked="" type="checkbox"/> CO2	1.860		11			100
<b>Background intensity:</b>		.000000	.000000	.000000	.000000	.000000
<b>Detector:</b>		Farad	Farad	Farad	Farad	Farad
<b>Ion Repeat:</b>		1	1	1	1	1

Preamp Autorange every 1 results  
\* enter 0 to autorange only once, in the beginning

Average every 1 results

☐ Fast Acquisition Mode

23. Analysis results will now be displayed in the "Analysis Results" page

Questor5 Client

Version: 3.30 Logout

Home Configuration Tuning Survey Scan **Analysis** Sequencer Results Event Log Compounds Help

Analysis Files [Analysis Method] Stream Evaluator Files Stream Evaluator Instrument Status

Test Analysis (Active) Valve: 1:No Valves

Results @ 08/31/2017 11:30:13

☒ Instrument ☒ Out Of Range

**Intensity Tags**

V1_I_18	0.016614152119
V1_I_28	0.009499258362 M
V1_I_32	0.039959009737
V1_I_40	0.009029956535
V1_I_44	0.004063971806

**Concentration Tags**

V1_C_H2O	26.35516630 %
V1_C_N2	11.48769034 %
V1_C_O2	51.74473813 %
V1_C_AR	7.63962352 %
V1_C_CO2	2.77277941 %

**Derived Tags**

**Analysis Results**